<u>Update instructions if using a pre-5.03 version of the Host PC</u> <u>Application with a Standalone Host PC (no Display PC connected)</u>


- 1) First, download the update zip file (e.g., "elcl_5.03.zip"). **Do not unzip it**. Save it on a USB drive. (the drive must use Fat32 format)
- 2) Then, on the Host PC, exit the Host PC Application by pressing Ctrl-Alt-Q on the Host PC keyboard. This will take you to the File Manager.
- 3) Insert the USB drive mentiond in step 1 to the Host PC.
- 4) On the Host PC, from the File Manager, click on the Configurations button:




5) Then, click on the System Update button:




6) Then, in the System Update dialog, click on the Choose File button:

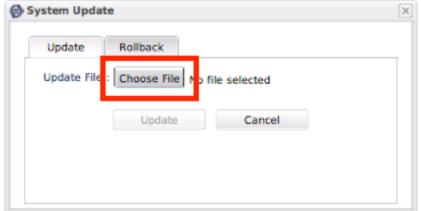


7) Then, in the Open dialog, click on the top bar marked "Look in:" and select the /fs/usb0 – this will be the USB memory stick that you have inserted into the Host PC:

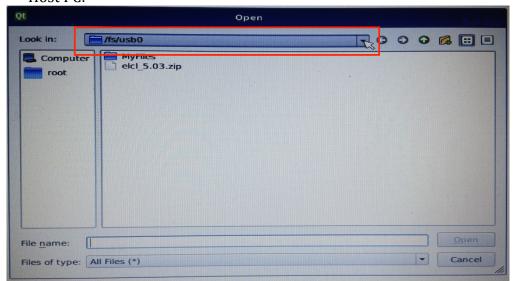

8) Select the update zip file and choose Open:

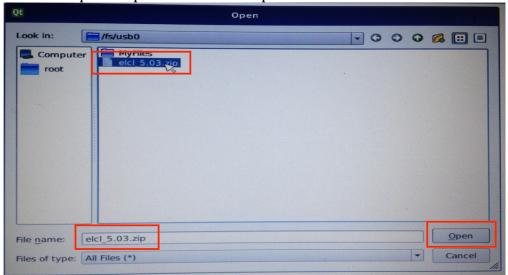

9) After the file has been selected, choose Update:

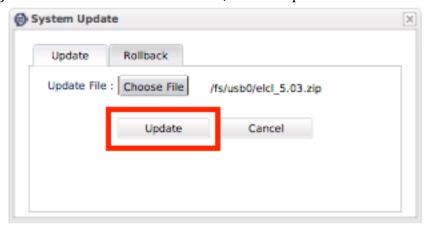

Wait until the host software is updated - please be patient as this process may take a few minutes to complete.

10) You can then select the /fs/usb0 device in the tree view of the File Manager and click the Eject button:

